Kalteinbettmittel

VariKleer

Best. Nr. 20-3590 (1 Satz)
Best. Nr. 20-3591 (1000 g Pulver)
Best. Nr. 20-3592 (500 ml Härter)

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Description:

VariKleer is an acrylic for general use that hardens crystal clear when cured in a pressure pot. If cured under normal pressure, the sample will become milky grey.

Instruction for use:

Specimens have to be clean and degreased prior to embedding.

☞ **Mix 2 Vol.-Parts of Powder with 1 Vol.-Part of Hardener Liquid in a small disposable container. Stir well and so that no air bubbles are stirred in.**

Processing time is about 3 min, curing time 10-15 min., depending on environmental temperature. In order to achieve crystal clear samples, VariKleer has to be cured in a pressure pot.

Max. curing temperature: 100°C

VariKleer can be used with every kind of moulds. In order to achieve crystal clear samples Silicon Rubber Mounting Moulds should NOT be used.

Caution:

Contains: Methyl methacrylate, Highly flammable • Irritant

R37/38: Irritating to respiratory system and skin. R43: May cause sensitisation by skin contact.

S16: Keep away from sources of ignition. No smoking. S23.3: Do not breathe vapor. S24: Avoid contact with skin.

S37: Wear suitable gloves. S9: Keep container in a well-ventilated place.